

# Courtesy Copy

SSD:TM:513

Rev. 8

Issue Date:  
11/13/2023

Implementation  
Date:  
11/27/2023

Page 1 of  
15

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

## Determination of Metals in Alcoholic Beverages by ICP-MS

---

### Scope and Application

This method will be used for the determination of metals in alcoholic beverages such as wines, distilled spirits, and malt beverages using inductively coupled plasma mass spectrometry (ICP-MS). This method is applicable to samples containing 0-100% alcohol by volume.

At the time of this revision, chromium, cobalt, nickel, copper, arsenic, selenium, cadmium, antimony, cesium, thallium, lead, and uranium are included.

Samples with high levels of solids, such as crème liquors, should not be analyzed using this method. A special sample preparation scheme is needed for samples of this type.

### Regulatory Tolerances

The tolerance levels of these elements are determined on the basis of regulations enforced by the Environmental Protection Agency (EPA), United States Food and Drug Administration (USFDA) and Alcohol and Tobacco Tax and Trade Bureau (TTB).

# Courtesy Copy

SSD:TM:513

Rev. 8

Issue Date:  
11/13/2023Implementation  
Date:  
11/27/2023Page 2 of  
15

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

## Levels and Limitations

### a. Instrument Levels and Limitations

<b>Instrument Levels and Limitations</b>			
	<b>Linear Range</b>	<b>Limit of Detection</b>	<b>Limit of Quantitation</b>
Cr <sub>52</sub>	0.5 to 50 ppb	0.02 ppb	0.07 ppb
Co <sub>59</sub>	0.5 to 50 ppb	0.01 ppb	0.04 ppb
Ni <sub>60</sub>	1.0 to 50 ppb	0.28 ppb	0.94 ppb
Cu <sub>63</sub>	1.0 to 50 ppb	0.19 ppb	0.64 ppb
Cu <sub>65</sub>	1.0 to 50 ppb	0.18 ppb	0.59 ppb
As <sub>75</sub>	0.5 to 50 ppb	0.01 ppb	0.05 ppb
Se <sub>82</sub>	0.5 to 50 ppb	0.02 ppb	0.08 ppb
Cd <sub>111</sub>	0.5 to 50 ppb	0.02 ppb	0.06 ppb
Sb <sub>121</sub>	0.5 to 50 ppb	0.04 ppb	0.14 ppb
Cs <sub>133</sub>	0.5 to 50 ppb	0.01 ppb	0.03 ppb
Tl <sub>205</sub>	0.5 to 50 ppb	0.01 ppb	0.02 ppb
Pb <sub>207</sub>	0.5 to 50 ppb	0.06 ppb	0.20 ppb
U <sub>238</sub>	0.5 to 50 ppb	0.01 ppb	0.03 ppb

# Courtesy Copy

SSD:TM:513

Rev. 8

Issue Date:  
11/13/2023Implementation  
Date:  
11/27/2023Page 3 of  
15

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

## b. Method Levels and Limitations

<b>Method Levels and Limitations *</b>			
	<b>Linear Range</b>	<b>Limit of Detection</b>	<b>Limit of Quantitation</b>
Cr <sub>52</sub>	10 to 1000 ppb	0.42 ppb	1.39 ppb
Co <sub>59</sub>	10 to 1000 ppb	0.22 ppb	0.71 ppb
Ni <sub>60</sub>	20 to 1000 ppb	2.85 ppb	9.41 ppb
Cu <sub>63</sub>	20 to 1000 ppb	1.95 ppb	6.42 ppb
Cu <sub>65</sub>	20 to 1000 ppb	1.79 ppb	5.89 ppb
As <sub>75</sub>	10 to 1000 ppb	0.28 ppb	0.92 ppb
Se <sub>82</sub>	10 to 1000 ppb	0.49 ppb	1.62 ppb
Cd <sub>111</sub>	10 to 1000 ppb	0.36 ppb	1.19 ppb
Sb <sub>121</sub>	10 to 1000 ppb	0.84 ppb	2.77 ppb
Cs <sub>133</sub>	10 to 1000 ppb	0.17 ppb	0.58 ppb
Tl <sub>205</sub>	10 to 1000 ppb	0.15 ppb	0.49 ppb
Pb <sub>207</sub>	10 to 1000 ppb	1.19 ppb	3.91 ppb
U <sub>238</sub>	10 to 1000 ppb	0.17 ppb	0.58 ppb
* Assuming a 1 to 20 dilution is performed			

# Courtesy Copy

SSD:TM:513

Rev. 8

Issue Date:  
11/13/2023Implementation  
Date:  
11/27/2023Page 4 of  
15

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

## Supplemental Documents

1. BAL:FORM:513-1 ICP-MS Daily Log
2. BAL:FORM:513-3 ICP-MS LCS Control Record
3. BAL:FORM:513-5 ICP-MS Daily Performance Record KED Mode (NLC\_4\_002)
4. BAL:FORM:513-6 ICP-MS Daily Maintenance (NLC\_4\_002)
5. BAL:FORM:513-7 Control Record for 5 ppb check
6. BAL:FORM:513-8 Thermo ICP-MS Semi Annual Maintenance (NLC\_4\_002)
7. BAL:FORM:302-4 ICP-MS Standards Log Record
8. NLC:FORM:514 ICP-MS Daily Maintenance Perkin Elmer Nexion (NLC\_4\_016)
9. TL:FORM:525-10 ICP-MS Daily Performance Record (NLC\_4\_016)
10. SSD:WG:513 Reagents and Sample Preparation and Handling for SSD:TM:513
11. SSD:WG:115 Procedure for Degassing Beer and Wine
12. NLC:WG:513-1 iCap-Q ICP-MS Instructions
13. NLC:WG:513-2 ICP-MS Semi-Annual Maintenance Schedule (NLC\_4\_002)

## Equipment

This method provides information for using the Perkin-Elmer NexION 2000 ICP-MS or the Thermo Scientific iCap Q ICP-MS

1. **Perkin-Elmer NexION 2000 ICP-MS** Inductively Coupled Plasma Mass Spectrometer or equivalent

ESI SC4 DXCi Autosampler or equivalent

### Instrument Conditions and Method Parameters

**Note:** Some of these values are approximations, some can change depending on the tuning of the instrument.

Nebulizer	Concentric PFA-ST or Glass with 400 $\mu$ L/min flow rate
Spray Chamber	Baffled Cyclonic
Sampler, Skimmer Cone and Hyper Skimmer Cone	Nickel or Platinum (Hyper Skimmer is aluminum)
RF Power	1600 W
Plasma gas flow rate	15.0 L/min
Nebulizer gas flow rate	0.85-1.1 mL/min
Auxiliary gas flow rate	1.0 – 1.20 L/min

# Courtesy Copy

SSD:TM:513

Rev. 8

Issue Date:  
11/13/2023Implementation  
Date:  
11/27/2023Page 5 of  
15

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

Sample uptake flow rate	1.0 mL/min
RPQ	0.25
Sweeps/reading	20
Readings/Replicate	1
Replicates	3
Dwell time	50 ms per AMU
Mode of analysis	Helium KED
Measurement Units	Counts per Second (CPS)

## 2. Thermo Scientific iCap Q ICP-MS Inductively Coupled Plasma Mass Spectrometer or equivalent

ESI SC4 DX Autosampler or equivalent

### Instrument Conditions and Method Parameters

**Note:** These are approximate values, some can change depending on the tuning of the instrument.

Nebulizer	Concentric PFA-ST or Glass with 400 $\mu$ L/min flow rate
Spray Chamber	Quartz Cyclonic spray chamber
Sample cone, Skimmer cone and Skimmer cone insert	Nickel or Platinum (Skimmer cone insert is nickel)
RF Power	1550 W
Plasma gas flow rate	14.0 L/min
Nebulizer gas flow rate	1.0 – 1.1 L/min
Auxiliary gas flow rate	0.80 L/min
Sample uptake	15 s
Spacing	0.2
Main runs	3

# Courtesy Copy

SSD:TM:513

Rev. 8

Issue Date:  
11/13/2023Implementation  
Date:  
11/27/2023Page 6 of  
15

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

Survey runs	1
Dwell time	0.03 s
Sweeps	10
Time per sweeps	0.63 s
Time per main run	6.3 s
Mode of analysis	KED
Measurement Units	Counts per Second (CPS)

**Signal Processing:** Applies only to Perkin Elmer NexION 2000 ICP-MS

Detector	Dual
Process Spectral Peak	Average
QID	On
Isotope Ratio Mode	Off
Blank Subtraction	After Internal Standard
Process Signal Profile	Average
Baseline Readings	0
Apply Smoothing Factor	5

## **Interferences**

1. Some metals, including arsenic, have strong Ar-Cl interferences that require correction compensation in the software.
2. Alcohol and other sources of carbon will interfere with metals analysis. Interference is minimized by preparation of standards in a similar carbon matrix. The addition of a small percentage of ethanol might be necessary to mitigate these carbon related interferences.
3. Some known interferences can be compensated by the use of correction equations in the processing method.

# Courtesy Copy

SSD:TM:513

Rev. 8

Issue Date:  
11/13/2023Implementation  
Date:  
11/27/2023Page 7 of  
15

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

## Elements Monitored

Element	Mass	Mode
Cr	52	KED
Co	59	KED
Ni	60	KED
Cu	63 65	KED KED
As	75	KED
Se	82	KED
Cd	111	KED
Sb	121	KED
Cs	133	KED
Tl	205	KED
Pb	207	KED
U	238	KED

## Internal Standards

The internal standard pairings at the time of this revision are:

Internal Standard	Elements grouped with this Internal Standard
Sc-45	Cr, Co, Ni, Cu (63, 65), As, Se
Y-89	Cd
In-115	Sb, Cs, Tl, Pb
Bi-209	U

## NOTE:

The internal standard pairings can be changed if necessary. Note any deviations.

<b>Courtesy Copy</b>	SSD:TM:513	Rev. 8
	Issue Date: 11/13/2023	Page 8 of 15
	Implementation Date: 11/27/2023	

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

## Reagents and Sample Preparation and Handling

Please refer to **SSD:WG:513** Reagents and Sample Preparation and Handling for SSD:TM:513.

## Procedures

The instrument should be optimized as per vendor instructions and/or internal guidelines.

## Example Sequence Tables

For the Thermo Scientific iCap Q ICP-MS, paired with the ESI SC4 DX Autosampler: Positions 1-21 on rack 1: use 50 mL conical tubes; on racks 2, 3 and 4 use 15 mL conical tubes. The Perkin-Elmer NexION 2000 ICP-MS is paired with the ESI SC4 DXCi Autosampler: Standard rack positions 1-9 use 125 mL bottles; Positions 101-121 on rack 1, Positions 201-260 on rack 2, positions 301-360 on rack 3 and positions 401-460 on rack 4 use 15 mL conical tubes.



# Courtesy Copy

SSD:TM:513

Rev. 8

Issue Date:  
11/13/2023

Implementation  
Date:  
11/27/2023

Page 9 of  
15

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

Example Sequence: This example is specific to the Thermo Scientific iCap Q ICP-MS, the sequence can be modified accordingly depending on the number of samples to be analyzed or any other situation.

Index	Label	Sample Type	Dilution Factor	Rack Number	Vial Numbers (Sample Position)	Survey Runs	Main Runs
1	Method Blank #1	UNKNOWN	1	1	7	1	3
2	Method Blank #2	UNKNOWN	1	1	7	1	3
3	Calibration Blank	Blk	1	1	1	1	3
4	0.5 ppb calibration Std	STD	1	1	2	1	3
5	1 ppb calibration Std	STD	1	1	3	1	3
6	5 ppb calibration Std	STD	1	1	4	1	3
7	25 ppb calibration Std	STD	1	1	5	1	3
8	50 ppb calibration Std	STD	1	1	6	1	3
9	Method Blank #3***	UNKNOWN	1	1	7	1	3
10	5 ppb Std Check (2nd Source) ***	QC	1	1	8	1	3
11	Method Blank #4	UNKNOWN	1	1	7	1	3
12	Unspiked LCS (unspiked Wine)*	QC	20	3	1	1	3
13	Method Blank #5	UNKNOWN	1	1	7	1	3
14	LCS-1* (spiked LCS)	QC	20	3	2	1	3
15	Method Blank #6	UNKNOWN	1	1	7	1	3
16	5 ppb Std Check 1 (2nd Source Accuracy)	QC	1	1	8	1	3
17	Method Blank #7	UNKNOWN	1	1	7	1	3
18	5 ppb Std Check 2 (2nd Source Precision)****	QC	1	1	9	1	3
19	Method Blank #8	UNKNOWN	1	1	7	1	3
20	5 ppb Std Check (2nd Source) ***	QC	1	1	8	1	3
21	Method Blank #9	UNKNOWN	1	1	7	1	3
22	Sample	UNKNOWN	20	3	3	1	3
23	Sample	UNKNOWN	40	3	4	1	3
24	Sample	UNKNOWN	50	3	5	1	3
25	Sample	UNKNOWN	100	3	6	1	3
26	Sample	UNKNOWN	200	3	7	1	3
27	Method Blank #10	UNKNOWN	1	1	7	1	3
n-1	5 ppb Std Check (2nd Source) ***	QC	1	1	8	1	3
n**	Method Blank #11	UNKNOWN	1	1	7	1	3

# Courtesy Copy

SSD:TM:513	Rev. 8
Issue Date: 11/13/2023	Page 10 of 15
Implementation Date: 11/27/2023	

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

Example Sequence: This example is specific to the Perkin-Elmer NexION 2000 ICP-MS, the sequence can be modified accordingly depending on the number of samples to be analyzed or any other situation. This is the view of the sequence after clicking “build run list.”

A/S Location	Sample ID	Measurement Action	Method	Sample Type
101	Unspiked Wine (Conditioning)*****	Run Diluted Sample	ssd tm 513_20x.mth	Sample
101	Unspiked Wine (Conditioning)*****	Run Diluted Sample	ssd tm 513_20x.mth	Sample
101	Unspiked Wine (Conditioning)*****	Run Diluted Sample	ssd tm 513_20x.mth	Sample
1	Method Blank #1	Run Sample	ssd tm 513_20x.mth	Sample
1	Method Blank #2	Run Sample	ssd tm 513_20x.mth	Sample
1	Method Blank	Run Blanks, Standards and Samples	ssd tm 513_20x.mth	Sample
1	Calibration Blank	Run Sample	ssd tm 513_20x.mth	Sample
9100	0.5 ppb Calibration Standard #1	Run Sample	ssd tm 513_20x.mth	Sample
9050	1.0 ppb Calibration Standard #2	Run Sample	ssd tm 513_20x.mth	Sample
9010	5.0 ppb Calibration Standard #3	Run Sample	ssd tm 513_20x.mth	Sample
9002	25.0 ppb Calibration Standard #4	Run Sample	ssd tm 513_20x.mth	Sample
9	50.0 ppb Calibration Standard #5	Run Sample	ssd tm 513_20x.mth	Sample
1	Method Blank #3***	Run Sample	ssd tm 513_20x.mth	Sample
7010	5 ppb Std Check (2nd Source) ***	Run Sample	ssd tm 513_20x.mth	Sample
1	Method Blank #4***	Run Sample	ssd tm 513_20x.mth	Sample
301	Unspiked LCS (unspiked Wine)*	Run Diluted Sample	ssd tm 513_20x.mth	Sample
1	Method Blank #5	Run Sample	ssd tm 513_20x.mth	Sample
302	LCS-1* (spiked LCS)	Run Diluted Sample	ssd tm 513_20x.mth	Sample
1	Method Blank #6	Run Sample	ssd tm 513_20x.mth	Sample
7010	5 ppb Std Check 1 (2nd Source Accuracy)	Run Sample	ssd tm 513_20x.mth	Sample
1	Method Blank #7	Run Sample	ssd tm 513_20x.mth	Sample
8010	5 ppb Std Check 2 (2nd Source Precision)****	Run Sample	ssd tm 513_20x.mth	Sample
1	Method Blank #8	Run Sample	ssd tm 513_20x.mth	Sample
7010	5 ppb Std Check (2nd Source) ***	Run Sample	ssd tm 513_20x.mth	Sample
1	Method Blank #9	Run Sample	ssd tm 513_20x.mth	Sample
303	Sample	Run Diluted Sample	ssd tm 513_20x.mth	Sample
304	Sample	Run Diluted Sample	ssd tm 513_40x.mth	Sample
305	Sample	Run Diluted Sample	ssd tm 513_50x.mth	Sample
306	Sample	Run Diluted Sample	ssd tm 513_100x.mth	Sample
1	Method Blank #10	Run Sample	ssd tm 513_20x.mth	Sample
7010	5 ppb Std Check (2nd Source) ***	Run Sample	ssd tm 513_20x.mth	Sample
1	Method Blank #11	Run Sample	ssd tm 513_20x.mth	Sample

# Courtesy Copy

SSD:TM:513

Rev. 8

Issue Date:  
11/13/2023

Implementation  
Date:  
11/27/2023

Page 11 of  
15

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

\* The unspiked LCS (unspiked Wine) will be placed before LCS 1 (spiked LCS). The spiked LCS does not have to be placed in any particular spot.

\*\* n: number of sequence lines/number of injections in sequence

\*\*\* Method Blank/2<sup>nd</sup> Source Check Standard can be placed in additional positions if necessary.

\*\*\*\* The 2<sup>nd</sup> source check for precision can be placed in any other position in the sequence.

\*\*\*\*\* The Unspiked wine (Conditioning) injections are placed at the beginning of the run only for the Nexion 2000 ICP-MS (Recommended by ESI).

## Quality Control

1. For the Thermo iCapQ ICP-MS every run should start with 1-2 method blanks to condition the system. When a run is stopped and re-started, it is best to condition the system with a method blank. These blanks are used to ensure all air bubbles have been removed from the pump tubing from the autosampler to the nebulizer. The method blank is the “blank” analyzed as a sample. The calibration curve begins with a calibration blank, which is subtracted from all samples and standards.
2. For the Perkin Elmer Nexion 2000 ICP-MS the sequence should start with 2-3 injections of wine to condition the system and cones. When a run is stopped and re-started, it is best to condition the system with a wine. These are used to ensure all air bubbles have been removed from the pump tubing from the autosampler to the nebulizer. The method blank is the “blank” analyzed as a sample. The calibration curve begins with a calibration blank, which is subtracted from all samples and standards.
3. A linear calibration curve will be used. The correlation coefficients of the calibration curves should be 0.99 or above for all elements.
4. The 5 ppb second source standard is analyzed as a check for accuracy. It will be analyzed in duplicate as a check for precision. The values should fall within  $\pm 30\%$  of the expected value (between 3 and 7 ppb) for the four monitored elements- Copper, Arsenic, Cesium and Lead.
5. The unspiked LCS (unspiked wine) and LCS (spiked LCS) will be analyzed, and the percent recovery will be checked and should be within 70 -130 % recovery for the four monitored elements Copper, Arsenic, Cesium and Lead.

# Courtesy Copy

SSD:TM:513

Rev. 8

Issue Date:  
11/13/2023

Implementation  
Date:  
11/27/2023

Page 12 of  
15

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

6. Run the 2<sup>nd</sup> source check standard after the calibration is completed, at least once every 10 samples and at the end of the sequence. Values for this standard should be within  $\pm 30\%$  of the expected value (between 3 and 7 ppb for all elements).
  
7. **Each occurrence of a failure for accuracy requires a CAR be initiated.** Refer to **SSD:QPD:202** for additional instructions.
  - a. If any of the four monitored elements in the 5 ppb second source standard for accuracy check fail then:
    - i. Re-run the 5 ppb second source standard. If all the values pass the second time, report all the results from the 5 ppb second source standard re-run and the sample data.
    - ii. If any value(s) fail a second time, do not report any results and pause the sequence (if possible). Inspect the different components of the instrument to see if a cause of the failure can be identified. Prepare a new 5 ppb second source standard or a 50 ppb multi-element second source stock standard and inject.

**Note:** If time allows and all QC parameters were met, the new 5 ppb second source or the 50 ppb multi-element second source stock standard preparation can be injected in the same sequence if the instrument was paused before injecting any samples. If time doesn't allow, run the new 5 ppb second source standard or 50 ppb multi-element second source stock standard in a new sequence. Also if using the Perkin Elmer Nexion 2000 ICP-MS the 5 ppb second source standard will be prepared from a 50 ppb multi-element second source stock standard by the ESI System.
    - iii. If all the new values pass then report all the results from the 5 ppb second source standard and report the sample data. If any of the values fail, do not report any results. Verify the instrument before shutdown, if no obvious reason for the failure can be identified shut down the instrument and inspect the different components to see if a cause of the failure can be identified.
    - iv. Refer to NLC-WG-513-1 for further trouble shooting (if using the Thermo Scientific iCap Q ICP-MS). Tag out the instrument and contact the co-principal analyst if the cause of the failure cannot be identified at the moment.

# Courtesy Copy

SSD:TM:513

Rev. 8

Issue Date:  
11/13/2023

Implementation  
Date:  
11/27/2023

Page 13 of  
15

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

- v. For precision failures the CAR can be logged in as an incident report (refer to SSD:QPD:310).
  - b. For 5 ppb second source check:
    - i. Values for this standard should be within  $\pm 30\%$  of the expected value (between 3 and 7 ppb for all elements).
    - ii. If any value is outside this range re-inject the second source, if all values are ok continue the analysis. If any value is still outside the allowed range pause the sequence (if possible). Inspect the different components of the instrument to see if a cause of the failure can be identified. Prepare a new 5 ppb second source standard or a 50 ppb multi-element second source stock standard (If using the Perkin Elmer Nexion 2000 ICP-MS) and inject. If all values are ok continue the analysis. If any value is still outside the allowed range shut down the instrument and inspect the different components to see if a cause of the failure can be identified.
    - iii. Refer to NLC-WG-513-1 (if using the Thermo Scientific iCap Q ICP-MS) for further trouble shooting. Tag out the instrument and contact the co-principal analyst if the cause of the failure cannot be identified at the moment.
8. It is recommended that the internal standard variation (ion counts) be  $100\% \pm 30\%$ . However, the results can be reported if the LCS recovery is within 70 – 130% and 2<sup>nd</sup> source standard checks are within  $\pm 30\%$  of the expected value.
9. A Method Blank should be run at least every 10 samples to check for carryover.
10. The following items should be submitted with sample data for the NexION 2000 ICP-MS.
  - a. Sequence template (if available)
  - b. Calibration curves and standards results
  - c. 5 ppb second source standard checks, calibration blank and method blanks
  - d. LCS and 5 ppb second source check data (accuracy and precision)

# Courtesy Copy

SSD:TM:513

Rev. 8

Issue Date:  
11/13/2023

Implementation  
Date:  
11/27/2023

Page 14 of  
15

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

11. If samples are analyzed with the Thermo Scientific iCap Q ICP-MS, the following items should be submitted with sample data:

- a. Instrument summary
- b. LabBook summary
- c. Method summary
- d. Sample list summary
- e. Calibration curves and standards results
- f. 5 ppb second source standard checks, calibration blank and method blanks
- g. LCS and 5 ppb second source check data (accuracy and precision)

## Sources of uncertainty:

Pipetting

Cross-contamination

## Calculations

$$LCS Recovery (\%) = \frac{Spiked LCS value (ppb) - Unspiked LCS value (ppb) \times 100}{Added amount (ppb)}$$

## Reporting Results

- Values should be reported to the nearest whole number (X ppb).
- Include a statement regarding the level the methodological measurement uncertainty is determined. For example, "The methodological measurement uncertainty is calculated on the basis of 100 ppb spike recoveries of the LCS elements, Cs, Cu<sub>63</sub>, Pb, and As."

## Safety Notes

Consult the MSDS for any chemicals used that are unfamiliar. All chemicals should be considered hazardous and avoid direct physical contact.

## Required Training, Certification, and Re-certification

1. In-house training by a certified chemist in the theory and operation of ICPMS, including software, maintenance, and troubleshooting.
2. Analyst has demonstrated competency after successfully obtaining 7 replicates of an unknown within the QC acceptance limits. (Prepare 7 spiked LCS samples and one unspiked LCS).
3. Chemists will be recertified periodically using proficiency test results and/or re-demonstration of competency.

# Courtesy Copy

SSD:TM:513

Rev. 8

Issue Date:  
11/13/2023

Implementation  
Date:  
11/27/2023

Page 15 of  
15

The colored ink stamp indicates this is a controlled document. Absence of color indicates this copy is not controlled and will not receive revision updates.

## Revision History

Rev. 1 – method converted from a “TMD” to a “TM”. Changes in bold italics represent edits from the draft method to the final TM. (4/11/2013)

Rev. 2 – included samples with higher alcohol content. Separated out the reagent and sample preparation and handling into a separate working guideline. Added “Related Documents” section.

Rev. 3 – changes to linear range, LOD, LOQ for copper (63 and 65) and nickel; addition of statement to narrative for measurement uncertainty in Reporting Results.

Rev. 4 – changes to Instrument linear range for Copper (63 and 65) and Nickel; addition to Quality Control Section regarding reporting the LCS.

Rev. 5 – Linear Ranges for Nickel and Copper (63 and 65) updated; Sources of Uncertainty section added.

Rev. 6 – Equipment section updated to reflect new instrumentation; Mode changed to KED in the Elements Monitored table; and updates made to Quality Control section.

Rev. 7 – Add additional supplemental documents. Add new Perkin Elmer Nexion 2000 ICP-MS, remove old Nexion ICP-MS. Edit instrument conditions and parameters for both instruments. Edit example sequence table and quality control sections.

Rev. 8 - Add additional supplemental documents. Edit instrument conditions, parameters, and signal processing for both instruments. Add information about carbon related interference mitigation. Edit example sequence tables and added clarifications to Quality Control section. Added information for competency requirements.